# AGAR, Bacteriological (7178)

# Intended Use

Agar, Bacteriological is a solidifying agent for use in preparing microbiological culture media.

# **Product Summary and Explanation**

Agar is a phycocolloid extracted from a group of red-purple algae, ususally *Gelidium* spp. Agar was first suggested for microbiological purposes in 1881 by Fannie Hesse.<sup>1,2</sup> By the early 1900's, agar became the gelling agent of choice over gelatin because agar remains firm at growth temperature for many pathogens and Agar is generally resistant to a breakdown by bacterial enzymes. The use of agar in microbiological media significantly contributed to the advance of microbiology, paving the way to study pure cultures.

Agar is a gel at room temperature, remaining firm at temperatures as high as 65°C.<sup>3</sup> Agar melts at approximately 85-91°C, a different temperature from solidification at 34-36°C. This property is known as hysteresis. Agar is generally resistant to shear forces; however, different agar may have different gel strengths or degrees of stiffness.

Specifications for Agar, Bacteriological include good clarity, controlled gelation temperature, controlled melting temperature, good diffusion characteristics, absence of toxic bacterial inhibitors, and relative absence of metabolically useful minerals and compounds. Agar, Bacteriological is recommended for clinical applications, auxotrophic studies, bacterial and yeast transformation studies, and bacterial molecular genetics applications.<sup>4,5</sup>

### **Principles of the Procedure**

Agar is typically used in a final concentration of 1 - 2% for solidifying culture media. Smaller quantities (0.05 - 0.5%) are used in media for motility studies (0.5% w/v), growth of anaerobes (0.1%) and microaerophiles.<sup>3</sup>

### **Precaution**

1. For Laboratory Use.

### **Quality Control Specifications**

Dehydrated Appearance: Powder is granular, homogeneous, free-flowing and creamy white beige.

Prepared Appearance (1.5% wt/vol): Prepared medium is very light amber to medium amber and slightly opalescent.

pH (2% Solution at 25°C): 6.0 - 7.5

**Gel Strength:** 760 - 850 g/cm<sup>2</sup>

**Expected Cultural Response:** Cultural response in Peptone Agar after incubation at 35°C for 18 - 24 hours incubation.

Microorganism	Response
Escherichia coli ATCC 25922	good to excellent growth
Staphylococcus aureus ATCC 25923	fair to good growth

### Test Procedure

Refer to appropriate references for specific procedures using Agar, Bacteriological.

# <u>Results</u>

Refer to appropriate references for test results.

### **Storage**

Store sealed bottle containing Agar, Bacteriological at 2 - 30°C. Once opened and recapped, place container in a low humidity environment at the same storage temperature. Protect from moisture and light by keeping container tightly closed.

### **Expiration**

Refer to expiration date stamped on container. Agar, Bacteriological should be discarded if not free flowing, or if the appearance has changed from the original color. Expiry applies to Agar, Bacteriological in its intact container when stored as directed.

### Packaging

Agar, Bacteriological	Code No.	7178A	500 g
		7178B	2 kg
		7178C	10 kg

#### References

- 1. Hesse, W. 1894. Uber die quantitative Bestimmung der in der Luft enthaltenen Mikroorganismen. Mit. a.d. Kaiserl. Gesh. Berlin. 2: 182-207.
- 2. Hitchens, A. P., and M. C. Leiking. 1939. The introduction of agar-agar into bacteriology. J. Bacteriol. 37:485-493.
- 3. Selby, H. H., and T. A. Selby. 1959. Agar. In Whister (ed.). Industrial gums, Academic Press Inc., New York, N. Y.
- 4. Sambrook, J., E. F. Fritsch, and T. Maniatis. 1989. Molecular cloning, a laboratory manual, 2<sup>nd</sup> ed. Cold Spring Harbor Laboratory Press, New York, N. Y.
- 5. Schiestl, R. H., and R. Daniel Geitz. 1989. High efficiency transformation of intact yeast cells using single stranded nucleic acids as a carrier. Current Genetics. 16:339-346.

### **Technical Information**

Contact Acumedia Manufacturers, Inc. for Technical Service or questions involving dehydrated culture media preparation or performance at (410)780-5120 or fax us at (410)780-5470.